



IKK β (I κ B Kinase β) Positive Control

Product Data Sheet

For Research Use Only, Not for use in diagnostic procedures

IKK β (I κ B Kinase β) Positive Control (Human, full length, recombinant enzyme expressed in Sf9.) Cat# CY-E1178-2

Lot No. 09C10

For 200 assays

(10 mUnits / μ L x 200 μ L)

Product Description: Human full length IKK β , containing a *N*-terminal GST tag and a C-terminal His tag, expressed in *Sf9*. Purified by GSH-agarose chromatography. The IKK β Positive control is designed to use for IKK α/β assay/Inhibitor Screening Kit [Cat# CY-1178]. The IKK β Positive Control should be added to the well at 10 m units/well. For instance, diluted positive control 1:10, use 10 μ L for 1 assay. Unused IKK β Positive control should be stored at -70°C.

Product Size: Full length IKK β : 2 units/200 μ L

Formulation: The IKK β Positive Control is supplied frozen in a buffer containing 20mM HEPES-KOH (pH 7.5), 1 % BSA, 2 mM DTT, 50mM NaCl, 0.03 % Brij35 and 50% glycerol.

Source: Human full length IKK β , containing *N*-terminal GST tag and a C-terminal His tag, expressed in *Sf9*.

Molecular Weight: IKK β demonstrates a single 105 kDa band by SDS-PAGE analysis.

Purity: IKK β is greater than 85% pure as determined by SDS-PAGE analysis.

Substrates: IKK β phosphorylates I κ B α at Ser32 and Ser36.

Inhibitors: BMS-345541, SC-514 and [5-(*p*-Fluorophenyl)-2-ureido]thiophene-3-carboxamide are known as a specific inhibitor for IKK β .

Unit Definitions: One unit is defined as the amount of kinase required to incorporate 1 nmol of phosphate into the I κ B α -S32/S36 peptide; KKKERLLDDRHDSGLDSMKDEEYE, per minute at 30°C.

Assay Conditions: Assay activity of IKK β in a 50 μ L reaction containing 20 mM Hepes KOH (pH 7.5), 5 mM MgCl₂, 1 mM DTT, 100 μ M [γ -³²P] ATP (1 μ Ci), and 4 μ g of I κ B α -S32/S36 peptide. Start the reaction by adding 10 μ L of the enzyme, diluted 50-fold in a buffer containing 20 mM Hepes KOH (pH 7.5), 1 mM DTT, 0.03 % Brij35. Incubate for 30 minutes at 30°C. The reaction is terminated by "spotting" a sample onto a phosphocellulose P81 filter paper disc, followed by washing extensively with 70 mM phosphoric acid to remove unincorporated radiolabel. Subsequently rinse the filters with ethanol and dry. The incorporated radioactivity on P81 filter is counted in a liquid scintillation counter

Storage and Stability: Stable for 12 months at -70°C from date of shipment. For maximum recovery of product, centrifuge the original vial after thawing and prior to removing the cap. Aliquot enzyme to avoid repeated freezing and thawing.

Related Products:

* IKK α and β Assay/Inhibitor Screening Kit: Cat# CY-1178



IKK β (I κ B Kinase β) Positive Control

Product Data Sheet

For Research Use Only, Not for use in diagnostic procedures

General References:

1. Ghosh, S., and Karin, M. (2002) *Cell* **109**, S81–S96
2. Schmitz, L. M., Bacher, S., and Kracht, M. (2001) *Trends Biochem. Sci.* **26**, 186–190
3. Naumann, M., and Scheidereit, C. (1994) *EMBO J.* **13**, 4597–4607
4. Brown, K., Gerstberger, S., Carlson, L., Franzoso, G., and Siabenlist, U. (1995) *Science* **267**, 1485–1488
5. Chen, Z., Haglen, J., Palombello, V. J., Melandri, F., Scherer, D., Ballard, D., and Maniatis, T. (1995) *Genes Dev.* **9**, 1586–1597
6. DiDonato, J., Mercurio, F., Rosette, C., Wi-Li, J., Suyang, H., Ghosh, S., and Karin, M. (1996) *Mol. Cell. Biol.* **16**, 1295–1304
7. Alkalay, I., Yaron, A., Hatzubai, A., Jung, S., Avraham, A., Gerlitz, O., Pashut-Levon, I., and Ben-Neriah, Y. (1995) *Mol. Cell. Biol.* **15**, 1294–1301
8. Karin, M., and Ben-Neriah, Y. (2000) *Annu. Rev. Immunol.* **18**, 621–663
9. DiDonato, J. A., Mercurio, F., and Karin, M. (1995) *Mol. Cell. Biol.* **15**, 1302–1311
10. Scherer, D. C., Brockman, J., Chen, Z., Maniatis, T., and Ballard, D. (1995) *Proc. Natl. Acad. Sci. U. S. A.* **92**, 11259–11263
11. Spencer, E., Jiang, J., and Chen, Z. J. (1999) *Genes Dev.* **13**, 284–294
12. Zandi, E., Rothwarf, D. M., Delhase, M., Hayakawa, M., and Karin, M. (1997) *Cell* **91**, 243–252
13. Hu, M. C., and Wang, Y. (1998) *Gene (Amst.)* **222**, 31–40
14. DiDonato, J. A., Hayakawa, M., Rothwarf, D. M., Zandi, E., and Karin, M. (1997) *Nature* **388**, 548–554
15. Mercurio, F., Zhu, H., Murray, B. W., Shevchenko, A., Bennet, B. L., and Rao, A. (1997) *Science* **278**, 860–866
16. Rothwarf, D. M., Zandi, E., Natoli, G., and Karin, M. (1998) *Nature* **395**, 297–300
17. Yamaoka, S., Courtois, G., Bessia, C., Whiteside, S. T., Weil, R., Agou, F., Kirk, H. E., Kay, R. J., and Israel, A. (1998) *Cell* **93**, 1231–1240
18. Li, Q., Van Antwerp, D., Mercurio, F., Lee, K. F., and Verma, I. M. (1999) *Science* **284**, 321–32535.
19. Li, Z. W., Chu, W., Hu, Y., Delhase, M., Deernick, T., Ellisman, M., Johnson, R., and Karin, M. (1999) *J. Exp. Med.* **189**, 1839–1845
20. Peters, R. T.; Liao, S.-M. (2000) Maniatis, T. *Mol. Cell* **5**, 513-522

PRODUCED BY

CycLex Co., Ltd.
1063-103 Terasawaoka
Ina, Nagano 396-0002
Japan
Fax: +81-265-76-7618
e-mail: info@cyclex.co.jp
URL: <http://www.cyclex.co.jp>

CycLex/CircuLex products are supplied for research use only. CycLex/CircuLex products and components thereof may not be resold, modified for resale, or used to manufacture commercial products without prior written approval from CycLex Co., Ltd.. To inquire about licensing for such commercial use, please contact us via email.